
YG

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Pressurization rotary drum filter

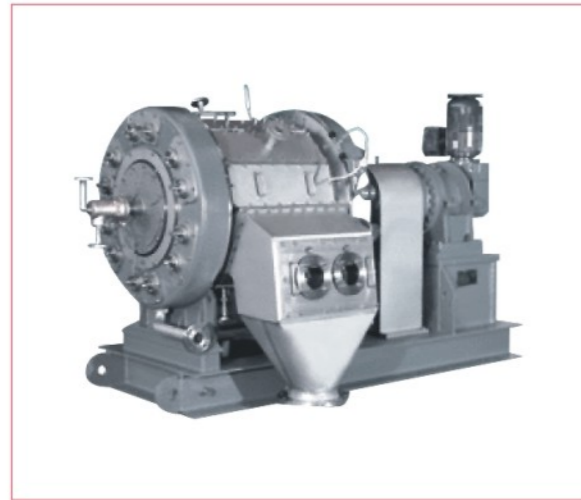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● Summary and application

The pressurization rotary drum filter, integrating washing, extracting, evaporating and drying, is a piece of closed equipment that realizes solid-liquid separation after continuously and automatically retaining the solid granules in the liquid and solid mixture by using porous rotary drum filtering media. It can be used to quickly carry out solid-liquid separation of various suspension materials with such features as continuous discharging, high handling capacity, good dehydrating effect, good operating environment, low floor space and high degree of automation. The product design conforms to the requirements in GMP and FDA.

● Technical Features

1. Washing cycle

The filter cakes with slightly extruded special structure and density can be washed in one or several phases. This kind of design in the rotary pressurization filter ensures that the minimum amount of washing media is used for replacement washing. The washing liquid is discharged through the liquid discharging pipe. The filter cakes are separated through the control head (depending on the washing phase selected) and return to the filter if necessary. The filter cakes go for drying and evaporating.

2. Filter cake discharging

Taking into consideration the operation of the full-automatic filter or the service life of the filter structure, it is very necessary to carry out simple and complete filter cake discharging. In the rotary pressurization filter, the filter cake will drop due to its own gravity. If the sticking force prevents the filter cake from falling, back blowing by air or gas or mechanical discharging by the scraper can serve as an aid to the discharging process. The filtering media are washed clean in the downstream washing chamber. The washing mechanism can use a pressure up to 25bar, which makes it possible to wash the filter clean from the retaining direction.

3. Continuous running

The previous pressurization filters almost adopt intermittent operation, but this machine adopts continuous operation, which conforms to modern requirements for continuous production. It displays its unique advantages because it can filter harmful and poisonous solids and solvents.

4. Perfect sealing performance

There is no need to worry about leakage even when it is used for handling filter pulps containing methyl and alcohol methylbenzene, or when it is necessary to adopt the volatile solvents as the detergent, or when materials of stinky smell are handled.

5. High efficiency

The rotary drum simultaneously completes the procedures such as filtering, filter cake washing, dehydrating and drying, discharging, etc. when it turns a circle.

6. Low water content of filter cakes

The water content of filter cakes is low because the compressed air, superheated steam or high temperature gas are used to dry the filter cakes in addition to the pressurization filter.

7. Safety

The potential risk of high speed rotary machinery will not occur because the rotation speed of the machine is low. Dangerous high-speed rotating machinery.

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
● 结构图 Block Diagram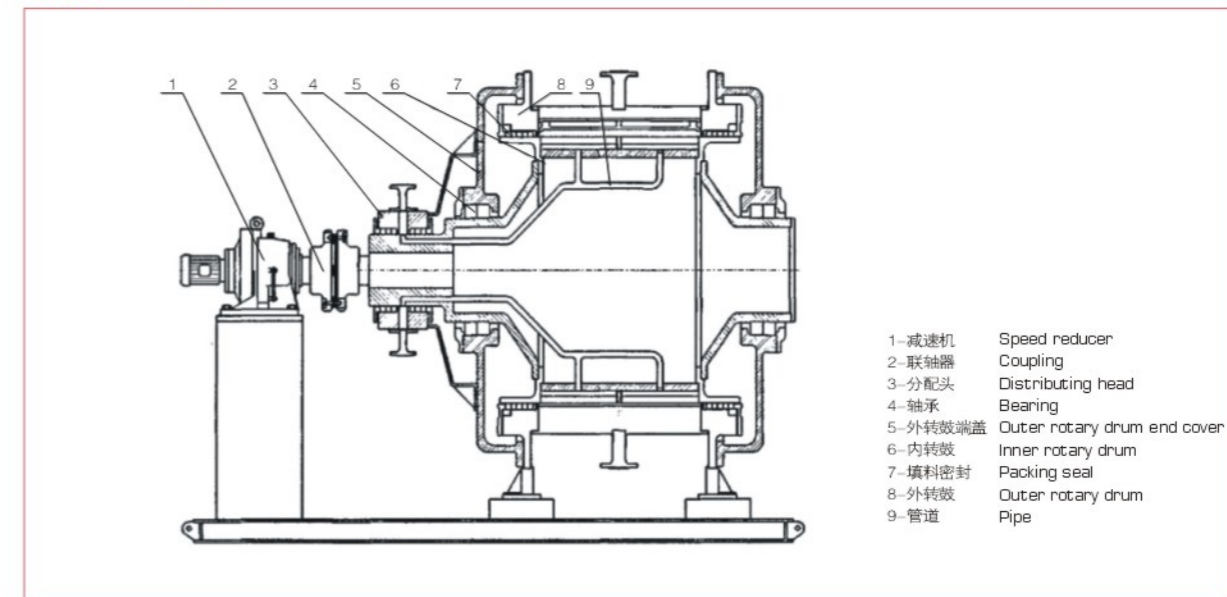

- 1-减速机 Speed reducer
- 2-联轴器 Coupling
- 3-分配头 Distributing head
- 4-轴承 Bearing
- 5-外转鼓端盖 Outer rotary drum end cover
- 6-内转鼓 Inner rotary drum
- 7-填料密封 Packing seal
- 8-外转鼓 Outer rotary drum
- 9-管道 Pipe